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


Single event spectroscopy on laser generated x-ray beams

C. Kern, M. Schnell, C. Spielmann, Institute of Optics and Quantum Electronics, University of Jena, Germany, (October 2010)



Introduction

The use of ultra intense laser pulses to excite plasma waves with a relativistic phase velocity is a possible route to the development of compact particle accelerators. Quasi-monoenergetic electron beams with energies from 0.1 to 1 GeV have been reliably generated [1,2]. In addition, these compact particle accelerators are sources of intense x-rays with peak brilliances comparable to "3rd generation" synchrotrons [4]. The electrons in the laser driven wakefield do not only experience a strong accelerating field but also a strong focusing field, forcing them to oscillate transversally in so called "betatron orbits" [3]. The pulse is perfectly synchronized with the driving laser pulse; the duration of the generated x-rays is of the same order of its pulse duration [5]. These unique properties offer the opportunity to investigate ultrafast processes in solids or warm dense matter using femtosecond x-ray diffraction or time-resolved absorption spectroscopy (NEXAFS / EXAFS).

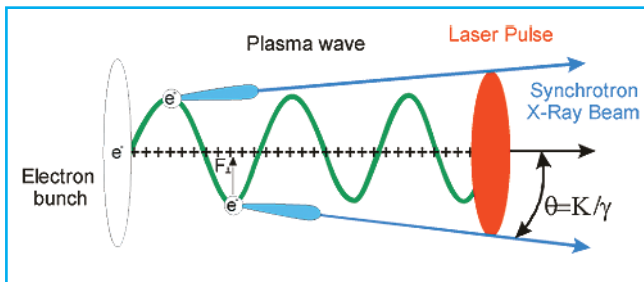


Fig. 1 Schematic of the production of short x-ray pulses by betatron oscillation inside a plasma

X-rays sources of this type are based on the betatron oscillation of a relativistic electron in a laser-produced plasma wave. Large numbers of these electrons are bunched together by the interaction with the driving field. Electrons inside each bunch are accelerated longitudinally with the propagating pulse. Additionally they experience the transverse electrostatic field within the wave, much as fast electrons experience periodic magnetic forces inside a synchrotron. The transverse acceleration leads to betatron oscillations. The result is a collimated beam of broadband radiation in the x-ray range which is emitted at femtosecond pulse lengths. The spectrum and maximum achievable photon energy is determined by the laser and gas parameters.

Application Note

Setup

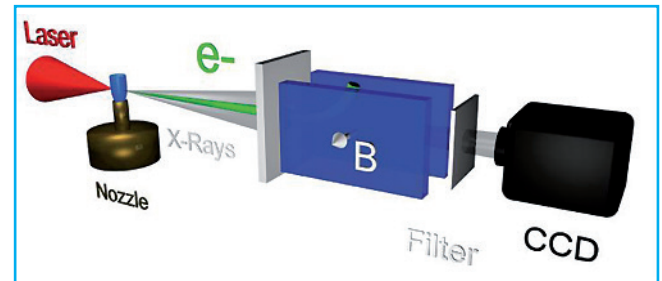


Fig. 2 Experimental setup for the detection and characterization of betatron radiation from laser-plasma-interaction in a He gas jet

As driving laser, we use the JETI laser system at the Institute of Optics and Quantum Electronics in Jena. This is a multistage CPA system, based on titanium doped sapphire, which emits at 800 nm central wavelength. For these experiments, pulse energies up to 0.7 J of 30 fs duration were used. This corresponds to peak intensities inside the target slightly above 10^{18} W/cm². The laser is focused into a supersonic Helium gas-jet (Fig. 2). The spectrum of the laser accelerated electrons was measured by deflecting them with a 0.5 T permanent magnet onto lanex screens (not shown in Fig. 2). The emitted x-rays are recorded by a CCD. This scheme allows to correlate the energy spectrum of the electrons with the energy of emitted photons for single shot experiments.

In order to characterize the emitted x-rays we use a large area back-illuminated CCD (iKon-L DO936N-MW-BN by Andor Technology). Thin filters were placed in front of the CCD to block radiation below 1 keV. The camera was calibrated for single photon event spectroscopy from well-defined classical x-ray sources. The count number of single events is assigned to a characteristic photon energy. The small pixel size ($13.5 \times 13.5 \mu\text{m}^2$) allows reasonable resolution, while the separation of neighboring events is still possible. The spatial resolution and high dynamic range can be useful when determining the spatial geometric properties of the source, e.g. by projecting the shadow image of a wire of known diameter onto the chip. In this way the source was determined to be smaller than $14 \mu\text{m}$ in diameter.

Single event spectroscopy on laser generated x-ray beams

C. Kern, M. Schnell, C. Spielmann, Institute of Optics and Quantum Electronics,
University of Jena, Germany, (October 2010)



Application Note

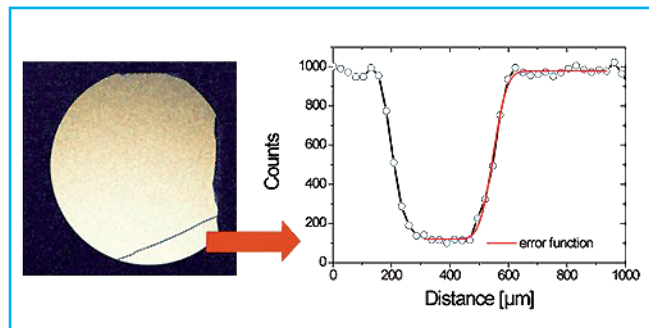


Fig. 3 Determination of the x-ray source size by evaluating the shadow thrown by a narrow wire projected on the CCD

Experimental results

So far, more than 10,000 single shot images were taken in this setup. From one image one can reconstruct a typical x-ray spectrum of the source. The results are shown in Fig. 4. We find photons of energies up to 10 keV, with a maximum in the spectrum around 2 keV. The low energy cutoff is given by the transmission of the used filters. Accounting for the filter function by deconvolution, the results agree reasonably well with simulations.

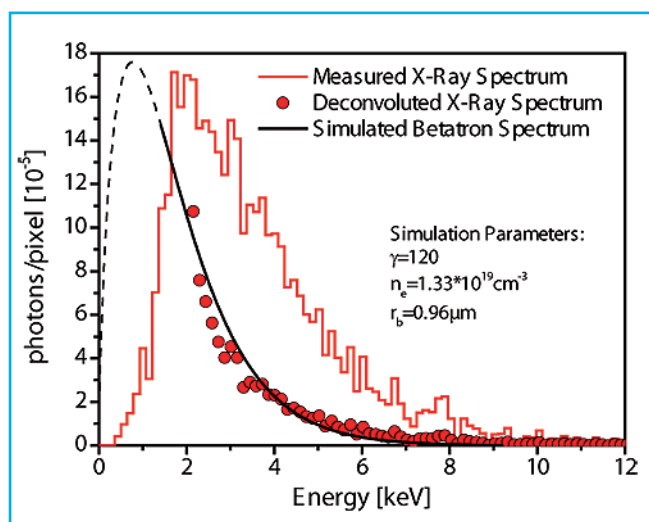


Fig. 4 X-ray spectrum of the betatron source

Summary & Outlook

First experiments have successfully demonstrated the generation of broadband betatron radiation with photon energies of several keV. In the future density shaped plasmas will be investigated to generate x-rays with even higher photon energies. The long term goal is the use of betatron radiation for ultrafast time-resolved absorption spectroscopy. This work is funded as a subproject of the SFB Transregio 18.

References

- [1] Leemans et al, Nat. Phys. 2 696 (2006)
- [2] Kneip et al., PRL 103, 035002 (2009)
- [3] Rousse et al., PRL 93 (2004)
- [4] Kneip et al., PRL 100, 105006 (2008)
- [5] K.Ta Phuoc et al., Phys. Plasmas, 14, 080701 (2007)

Contact

Michael Schnell
Institute of Optics and Quantum Electronics
University of Jena
Max-Wien-Platz 1
07743 Jena
Germany
phone: +49 (0)3641 947268
fax: +49 (0)3641 947232
email: michael.schnell.1@uni-jena.de
www: <http://www.physik.uni-jena.de/qe/start.html>